

Fire Monkey, a spin column HMW-DNA extraction kit

Dr. Georgios Patsos, CSO & kit inventor



Spin columns for long reads

Rapid & user-friendly extraction of HMW-DNA or RNA from mammalian cells or bacteria



- Spin columns NAIPs are fast & user-friendly
- High g-force applied guarantees extraction speed/efficiency
- However, high g-force breaks DNA
- Fire Monkey patented chemistry & matrix overcomes that
- Protocols remain fast and user-friendly
- Filtration removes small fragments-no need for size selection
- Spinning homogenizes long fragment distribution
- 100kb+ average strand length-no need for fragmentation
- Extract is library-ready
- Parallel RNA extraction possible

~1hr



Kit



Equipment

Sample type	Throughput (Gb)	N50 (kb)
Gram negative	25.63	45.8
Gram positive	15.7	50.3
White blood cells	21.1	56

LSK109/MinION at $\geq Q7$

From sample to DNA in **1h**

Spin column
Standard format

> 100kb
Average extract size



FIRE MONKEY
HMW-DNA
extraction kit

Bacteria cells throughput up to **~26 Gb**

10-20kb
Size selection Cut-off

N50 mammalian cells up to **62 kb**

Automation

Library ready DNA for multiplexed long reads


Automation platform






Fluent 780 with integrated centrifuge and heated shakers

- Tecan developed an automated platform for HMW-DNA extraction utilising Fire Monkey chemistry & matrix
- Plate centrifuge proof-of-concept
- Bacteria were processed using Fire Monkey 96-well filter plates vs Fire Monkey columns
- Generated extracts with very similar extraction yield and average strand length
- Fluent 780 installation and full workflow optimization is next

Filter plate proof of concept

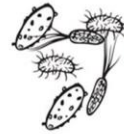

0.5 & 1 billion *E coli*

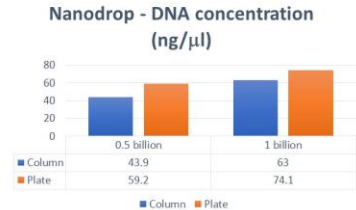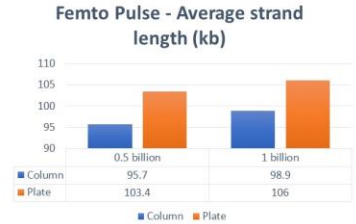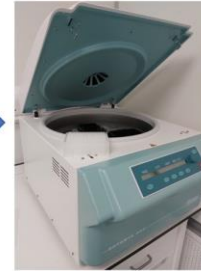

Automation/multiplexing applications



Updated extraction protocols



Efficient resuspension > major DNA recovery boost > more throughput

Needle cell resuspension-FM version 8

New protocol benefits




~1hr

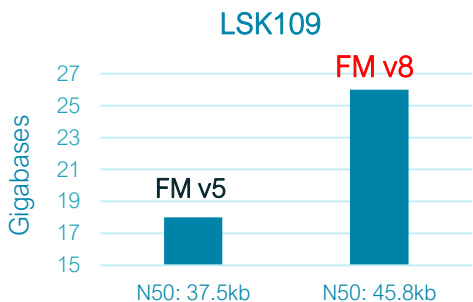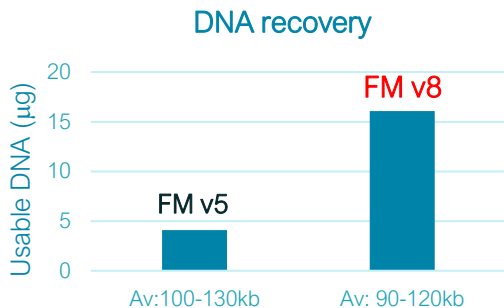

- Needle cell resuspension prior to Proteinase K lysis boosts DNA recovery
- New protocol version (v8) extracts ~4x more usable DNA than previous version (v5)
- DNA not 'too short/not too long' generates maximum throughput at a high N50
- Critical starting concentration (100ng/µl) achieved
- More DNA=more runs, or more loads per library
- More homogeneous HMW-DNA lowers \$/Gb price

N: NEB
C: Circulomics
R: RevoluGen
Q: Qiagen
P: Promega
Z: Zymo Research

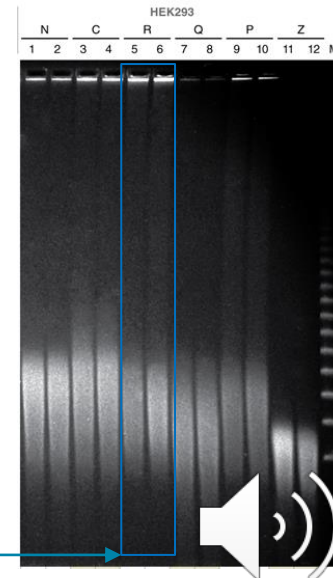

Old FM v5 version published by NEB

Thank you

The content in this presentation should not be reproduced without permission of the speaker. Oxford Nanopore Technologies, the Wheel icon, EPI2ME, Flongle, GridION, Metrichor, MinION, MinKNOW, Plongle, PromethION, SmidgION, Ubik, and VolTRAX are registered trademarks of Oxford Nanopore Technologies in various countries. All other brands and names are the property of their respective owners.

© 2020 Oxford Nanopore Technologies. All rights reserved.
Oxford Nanopore Technologies products are currently for research use only.

